



Smart Stir

Manual



Model


SS-100

List of Parts

Item	Picture	Description
Smart Stir Base	 A white, rounded rectangular base with a central opening and a small blue circular button on the front.	Holder Size : 1.06 inches (27 mm) in Diameter
Stirring Bar	 A white, cylindrical magnetic stirring bar.	<ul style="list-style-type: none">• Magnetic Stirring Bar for use with Smart Stir• Size : 0.12 x 0.28 inch (3 x 7 mm)
Iron Wire	 A thin, silver-colored metal wire.	For taking out the stirring bar from the vial
Micro USB Charging Cable	 A white charging cable with a standard USB-A connector on one end and a Micro USB-B connector on the other.	For charging Smart Stir

Operation

- 1 Fill the vial with your water sample and add the reagent according to the instructions of the test kit manufacturer.
- 2 Put the magnetic stirring bar in the vial.
- 3 Place the vial in the Smart Stir Base and press the power button.
- 4 Press the power button again to stop stirring. Or it will stop automatically after 5 minutes.
- 5 Don't forget to remove the stirring bar with the iron wire before emptying the vial.

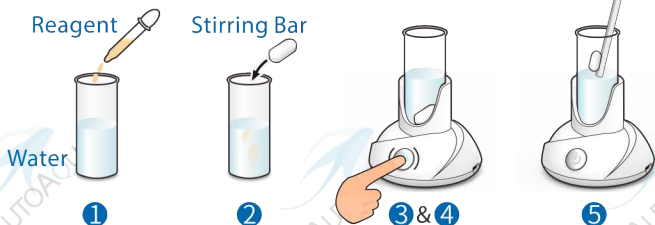

Notice

1. Please charge Smart Stir before first use.
2. Please fill the vial before starting stirring.
3. Please take out the stirring bar with the iron wire first before emptying the vial.
4. Smart Stir will stop operation automatically after five minutes.
5. The power button will flash blue when the battery is running low.
6. Please do not charge Smart Stir with a fast charger or it may be damaged.

Specification

- Charging : USB 5V
- Charging Time : 120 minutes
- Battery Life : 8 hours
- Smart Stir Base Size : 2.77 x 2.77 x 2.34 inches (70.5 x 70.5 x 59.5 mm)
- Holder Size : 1.06 inches (27 mm) in Diameter
- Stirring Bar Size : 0.12 x 0.28 inch (3 x 7 mm)

Limited Warranty

1. Limited warranty is provided by AUTOAQUA only if your AUTOAQUA Product ("Product") was purchased from an authorized distributor or an authorized reseller.
2. All warranty repairs or replacement are performed by the authorized distributor or authorized reseller in the region where the Product was first shipped by AUTOAQUA.
3. The limited warranty is valid for a period of 1 year under proper installation and normal use from the date of the original purchase listed on the proof of purchase or sales receipt.
4. In the event of a limited warranty claim, proof of purchase or sales receipt (issued to the original purchaser) is required.

Purchase Date

(Y/M/D)

Warranty Period

One year from the purchase date

Place of Purchase

(Store name, phone, email, address)

Service

AUTOAQUA Technologies Co., Ltd.
7F-7., NO.229, Fuxing 2nd Rd., Zhubei
City, Hsinchu County, Taiwan
E-mail : service@autoaqua.com.tw
www.autoaqua.com.tw

* Please go to AUTOAQUA website to get the distributor information.

